

Clinet® 303

Ready-to-use disinfecting spray



Effective immediately !

- Ready-to-spray
- Disinfects all surfaces
- No contact with the hands
= no irritation risks
- No rinsing

Clinet® 303 is always near at hand. Its solution enables highly effective and quick disinfection.

Ready-to-spray, **Clinet® 303** covers very large surfaces and reaches places not reachable by wiping such as wall angles, various asperities...

Clinet® 303 guarantees excellent laboratory hygiene.

Applications

Disinfection and cleaning of large surfaces:

- In laboratories
- In hospitals
- In offices

How to use

- Spray at about 30 cm from the surface to disinfect
- Dose : $\pm 40 \text{ g/m}^2$ (appears quite humid) for 1 L
- Wipe the surface
- Do not rinse

Composition

- Isopropyl alcohol
- Glutaric aldehyde
- Dimethyl-didecyl ammonium chloride

Precautions

- Keep out of reach of children
- Do not swallow
- Avoid contact with the eyes (irritant) and the skin
- Inflammable
- Not for food contact surfaces



For more info, the safety data sheet is available on request.

Active on

Bactericidal: NF T 72.151- NF T 72.171

Antibacterial: NF T 72.190

Technical specifications

Clinet® 303		
Reference	310 303	
Dispenser	Spray bottle 1 L	
Box	Box of	6
	Size	28 x 19 x 27.5 cm
	Gross weight	6.3 kg

Clinet® Range



Clinet® 101

Disinfecting wipes

Ref.: 310 101



Clinet® 202

Disinfecting wipes
in individual pack

Ref.: 310 202




Clinet® Anti-Algae

Liquid for inhibiting
algae growth

Ref.: 310 400

Discover our complete range for microbiology

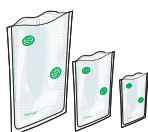

BagFilter®
Sterile filter bags



DiluFlow®
Gravimetric dilutors



BagMixer®
Lab blenders



FlexiPump®
Peristaltic pumps



easySpiral Dilute®
Automatic diluter and platers



Scan®
Automatic colony counters
Inhibition zone readers

ambifood®

Rua Dominguez Alvarez, 44, 4.16
4150-801 Porto Portugal

TEL: +351 229 962 069

FAX: +351 229 940 082

geral@ambifood.com

www.ambifood.com