

Clinet® Anti-Algae

Liquid for inhibiting algae growth



Efficiency for lab baths

- For water baths hygiene
- Reduces the water consumption
- Change the water every month only
- Easy-to-use thanks to its designed bottle

Clinet® Anti-Algae maintains and cleans the water baths and ultrasonic trays.

Applications

Prevents algae for water baths, pool baths...

How to use

- Use it diluted at 0.1 %:
1 dose of 20 mL for 20 L of demineralised water
- Renew the necessary dose every week
- Change the water bath every month

Composition

Ammonium quaternary, calcium, preservative

Precautions

- Use exclusively demineralized or deionized water because the algicide action is diminished by hard water.
- Do not mix with anion-active detergents and/or strong oxydants (such as bleach...)
- In case of contact with the eyes, rinse abundantly with water and consult a specialist
- Do not swallow
- Keep out of reach of children



For more info, the safety data sheet is available on request.

Technical specifications

Clinet® Anti-Algae	
Reference	310 400
Box of	1 L
Size	34 x 25.5 x 18.5 cm
Gross weight	1.1 kg

Clinet® range



Clinet® 101

Disinfecting wipes

Ref.: 310 101



Clinet® 202

Disinfecting wipes
in individual pack

Ref.: 310 202




Clinet® 303

Ready-to-use
disinfecting spray

Ref.: 310 303

Discover our complete range for microbiology

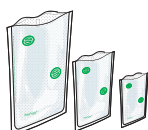

BagFilter®
Sterile filter bags



DiluFlow®
Gravimetric dilutors



BagMixer®
Lab blenders



FlexiPump®
Peristaltic pumps



easySpiral Dilute®
Automatic diluter and platers



Scan®
Automatic colony counters
Inhibition zone readers

ambifood®

Rua Dominguez Alvarez, 44, 4.16
4150-801 Porto Portugal

TEL: +351 229 962 069

FAX: +351 229 940 082

geral@ambifood.com

www.ambifood.com